## **New Product Highlights**

# 2,6-Difluoro-4-[2-(phenylsulfonylamino)ethylthio]phenoxyacetamide (PEPA): A novel, allosteric modulator of AMPA glutamate receptors Exclusively available from Sigma-RBI

The α-amino-3-hydroxy-5-methylisoxazole-4-propionic acid (AMPA) glutamate receptor subtype is a hetero-oligomer that is formed from combinations of four subunits referred to as GluR1-4. Each of these subunits exists in flip and flop isoforms that are produced by alternative mRNA splicing [1]. Two well-known and structurally distinct modulators of the AMPA receptor, **aniracetam** (Prod. No. <u>A 9950</u>) and **cyclothiazide** (Prod. No. <u>C 9847</u>) potentiate excitatory synaptic transmission by preferentially modulating the flop and flip splice variants respectively [2] and possess therapeutic potential as cognitive enhancers [3].

Sigma-RBI is pleased to introduce **2,6-difluoro-4-[2 phenylsulfonylamino)ethylthio]phenoxy-acetamide** (PEPA, Prod. No. **D 8941**), a structurally novel, selective, allosteric modulator of AMPA glutamate receptors. Using two-electrode voltage clamp recordings, PEPA potentiated glutamate-evoked currents in *Xenopus* oocytes by attenuating the extent of AMPA receptor desensitization expressing flop isoforms of GluR3 and GluR4 subunits, exhibiting an EC<sub>50</sub> value of approximately 50 μM. PEPA also reduced AMPA receptor desensitization in HEK293 cells transfected with all four GluR subunits, as measured by whole cell patch clamp recordings [3-5].

Although PEPA and aniracetam have similar subunit and splice-variant selectivity, at a concentration of  $10 \mu M$  PEPA was 100 times more potent than aniracetam in enhancing AMPA receptor currents in

Xenopus oocytes. PEPA and cyclothiazide have opposite splice-variant selectivity and at 25 μM PEPA potentiated glutamate responses in Xenopus oocytes to a much greater extent than cyclothiazide at the same concentration [3]. PEPA (1, 3, 10 mg/kg/day for 10 days) also ameliorated ischemia-induced impairment of performance produced by occlusion of the middle cerebral artery in rats, as assessed by Morris water maze test. In contrast, comparable doses of aniracetam did not significantly improve performance [6].


PEPA is therefore a structurally novel allosteric modulator of AMPA glutamate receptors and appears to be a highly selective tool with which to investigate the physiology of GluR subunits and their involvement in cognitive processes.

Sold under exclusive license from Nippon Suisan Kaisha, Ltd.

#### References

- 1. Sommer, B., et al., Science, 249, 1580-1585 (1990).
- 2. Partin, K.M., et al., J. Neurosci., 16, 6634-6647 (1996).
- 3. Sekiguchi, M., et al., J. Neurosci., 17, 5760-5771 (1997).
- Sekiguchi, M., et al., Br. J. Pharmacol., 123, 1294-1303 (1998).
  Sekiguchi, M., et al., Br. J. Pharmacol., 136, 1033-1041 (2002).
- Sekiguchi, M., et al., B. J. Frialmacol., 136, 1033-1041 (20
  Sekiguchi, M., et al., Neuroreport, 12, 2947-2950 (2001).

### **New Sigma-Aldrich On-Line Explorers**

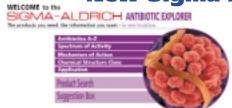

### **The Sigma-Aldrich Antibiotic Explorer**

Explore the industry's most comprehensive collection of Antibiotics with this new, easy-to-use web-based tool.

- Over 400 Antibiotics available on-line.
- Categorized according to Antibiotic applications.
- ◆ Monthly specials on the Antibiotics researchers use daily.
- Quick access to datasheets, certificates of analysis, material safety data sheets and technical service.

Visit the Antibiotic Explorer today at: <a href="mailto:sigma-aldrich.com/antibiotics">sigma-aldrich.com/antibiotics</a>



#### The Sigma-Aldrich Peptide Explorer

Looking for peptide products has been made easy.

- Products are conveniently categorized into 30 welldefined peptide families and sorted alphabetically.
- Search tips are provided to help you locate the peptide you need.
- Quick access to datasheets, certificates of analysis, material safety datasheets and technical service.

Visit the Peptide Explorer today at: sigma-aldrich.com/peptides

