

Product Information

Deoxyribonuclease I from bovine pancreas

Sigma Type II-S, lyophilized powder, protein $\geq 80\%$, ≥ 2000 units/mg protein**D4513**

Product Description

CAS Registry Number: 9003-98-9

Enzyme Commission (EC) Number: 3.1.21.1

Synonyms: DNase I, Deoxyribonuclease A, Deoxyribonuclease 5'-oligonucleotidohydrolase

Deoxyribonuclease I (DNase I) is an endonuclease that cleaves DNA by preferentially acting on phosphodiester bonds adjacent to pyrimidines, to produce polynucleotides with terminal 5'-phosphates. A tetranucleotide is the smallest average digestion product. In the presence of Mg^{2+} ions, DNase I attacks each strand of DNA independently and the cleavage sites are random. If Mn^{2+} ions are present, both DNA strands are cleaved at approximately the same site.¹ DNase I hydrolyzes single-stranded DNA, double-stranded DNA, and chromatin (the reaction rate is restricted by DNA association with histones).

DNase I is found in most cells and tissues. In mammals, the pancreas is one of the best sources for the enzyme. Pancreatic DNase I was the first DNase to be isolated. The calculated molecular mass is 30,072 Da. DNase I exists as a mixture of glycoproteins with two disulfide bridges.²

Bovine pancreatic DNase I contains four chromatographically distinguishable components, labeled A, B, C, and D.³ The molar ratios of A:B:C in a pancreatic extract are 4:1:1. Only minor amounts of D are found. Forms A and B differ in carbohydrate content (see Table below).⁴

Carbohydrate Content⁴

Carbohydrate / Form	A	B	C
N-Acetylglucosamine	2	3	2
Mannose	6	5	5
Sialic Acid	-	1	-
Galactose	-	1	-

Form C differs from Forms A and B by having one less His and one more Pro, and in the carbohydrate chain.⁴

DNase I is used to remove DNA from protein and nucleic acid samples, and to nick DNA as a first step to incorporate labeled bases into DNA. Several theses⁵⁻¹⁰ and dissertations¹¹⁻²⁶ have cited use of product D4513 in their protocols.

Isoelectric points:²

- A: 5.22
- B: 4.96
- C: 5.06
- D: 4.78

Optimal pH: 7-8

Extinction Coefficient: $E_{280}^{1\%} = 11.1$

Activators

- DNase I has an absolute requirement for divalent metal cations.
- The most commonly used divalent metal cation is Mg^{2+} .^{27,28}
- However, Mn^{2+} , Ca^{2+} , Co^{2+} , and Zn^{2+} will activate DNase I.²⁷⁻²⁹
- 5 mM Ca^{+2} will stabilize DNase I against proteolytic digestion.³⁰
- 0.1 mM Ca^{+2} is needed to reduce the rate of inactivation by one-half.³⁰

Inhibitors

There is no general inhibitor specific for DNase I.^{27,28} Citrate inhibits Mg^{2+} -activated DNase I, but not Mn^{2+} -activated DNase I.

- 2-Mercaptoethanol (the reduced enzyme is inactive, but can be reactivated in the presence of Ca^{2+} or Mg^{2+} ions)²⁸
- Chelators (such as EDTA, EGTA)
- Sodium dodecyl sulfate (SDS)³⁰
- Actin³²

MERCK

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Product

This product is chromatographically purified from bovine pancreas. The purification procedure is not selective for any form (A, B, C, or D) of DNase I. D4513 is supplied as a lyophilized powder, containing CaCl_2 . This product is 0.2 μm sterile-filtered. Each vial contains ~11 mg of protein.

Protein content: $\geq 80\%$ (Biuret)

Specific activity: ≥ 2000 Kunitz units/mg protein

Unit Definition:³³

- One Kunitz unit will produce a ΔA_{260} of 0.001 per minute per mL at pH 5.0 at 25 °C, using DNA, Type I or III, as substrate, with $[\text{Mg}^{2+}] = 4.2 \text{ mM}$.
- This enzyme assay reaction is performed in 95 mM acetate buffer, pH 5.0, at 25 °C, containing 4.75 mM Mg^{2+} and 1.9 mM Ca^{2+} , in a 3 mL reaction.

Contaminants:

- Protease: $\leq 0.005\%$
- RNase: $\leq 0.01\%$
- Chymotrypsin: $\leq 0.01\%$
- Endotoxin-tested

Preparation Instructions

This enzyme is soluble in 0.15 M NaCl at 5 mg/mL.

Storage/Stability

DNase I retains activity for at least three years, when unopened and stored long-term at the recommended temperature, -20°C .

Solutions of DNase I at 10 mg/mL in 0.15 M NaCl may lose $<10\%$ of its activity when stored for a week in aliquots at -20°C . The same solutions stored in aliquots at $2-8^\circ\text{C}$ can lose $\sim 20\%$ activity.

DNase I remains active in solution between pH 5-7 up to 60°C for at least five hours. A 1 mg/mL solution in acetate buffer (pH 5.0) or Tris buffer (pH 7.2) loses activity at the rate of 6% per hour. At 68°C , DNase I loses activity in <10 minutes.

References

1. Sambrook, J. *et al.*, *Molecular Cloning: A Laboratory Manual*, 2nd edition. Cold Spring Harbor Laboratory Press (Cold Spring Harbor, NY), Vol. 2, p. 5.83 (1989).
2. Liao, T.-H. *et al.*, *J. Biol. Chem.*, **248(4)**, 1489-1495 (1973).
3. Kim, H.S., and Liao, T.-H., *Anal. Biochem.*, **119(1)**, 96-101 (1982).
4. Salnikow, J. *et al.*, *J. Biol. Chem.*, **245(21)**, 5685-5690 (1970).
5. Ratanshi, Imran, "Evaluating the Capacity of Adult Human Mesenchymal Stem Cells to Differentiate into Insulin-Producing Cells *In Vitro*". University of Calgary, M.Sc. thesis, p. 82 (2008).
6. Govinder, Lerisa, "Development of novel T cell assays and assessment of immune recognition to latency associated *M. tuberculosis*-specific antigens Rv2660 and Rv2659". University of Cape Town, M.Sci. thesis, p. 90 (2010).
7. Chung, Cheong Kwet Choy Kwong, "Characterisation of the T cell responses induced by BCG in infants over the first year of life". University of Cape Town, M.Sci.Med. thesis, p. 109 (2011).
8. Garg, Aman, "Computational Agent-Based Modeling of Surgical Vocal Fold Injury and Repair". McGill University, M.Sc. thesis, p. 15 (2018).
9. Hündler, Manuel, "Isolating the iron storage protein ferritin from human brain tissue and characterizing its iron core via electron microscopic techniques". Graz University of Technology, M.Sc. thesis, p. 71 (2019).
10. Mahendralingam, Mathepan, "Lineage-restricted metabolic identities of mammary epithelial cells". University of Toronto, M.Sc. thesis, pp. 22, 28 (2019).
11. Hoppstädter, Jessica, "Differential expression of pro- and anti-inflammatory mediators in pulmonary macrophages upon Toll-like receptor activation". Universität des Saarlandes, Dr. rer. nat. dissertation, p. 111 (2010).
12. Laranjeira, Cátia Susana Torres, "*In vivo* Identification of Neural Stem Cells in the Enteric Nervous System". University College London, Ph.D. dissertation, p. 105 (2010).
13. Copley, Michael Rebin, "Regulation of Developmental Changes in Hematopoietic Stem Cell Self-Renewal". University of British Columbia, Ph.D. dissertation, p. 39 (2013).

14. Cheng, Chieh-Yang, "Understanding Prostate Cancer Pathogenesis". Cornell University, Ph.D. dissertation, pp. 59, 147 (2014).
15. Mansour, Tamer Ahmed, "Determining the Role of Irf6 in T Cell Development and Functional Commitment". Michigan State University, Ph.D. dissertation, p. 99 (2014)
16. Noges, Laura Evelyn, "Clarification of the Role of Host DNA as a Mediator of Vaccine Adjuvant Activity". University of Colorado, Ph.D. dissertation, p. 33 (2015).
17. Cosgrove, Megan Elizabeth, "The Role of JAK2^{V617F}-positive Endothelial Cells in Aberrant Haemostasis and Thrombosis". University of York, Ph.D. dissertation, p. 87 (2016).
18. Roberts, Allison Whitney, "Resident Macrophages are Locally Programmed for Silent Clearance of Apoptotic Cells". University of California Berkeley, Ph.D. dissertation, p. 20 (2016).
19. Timošenko, Elina, "Tryptophan catabolism and amino acid transporter reprogramming in the tumour microenvironment". University of Oxford, Ph.D. dissertation, p. 93 (2016).
20. Bilbao, Begoña Cánovas, "Role of p38 MAPK in breast cancer". Universitat de Barcelona, Ph.D. dissertation, p. 62 (2017).
21. Hume, Robert David, "Recapitulating mammary gland development and breast cancer cell migration in vitro using 3D engineered scaffolds". University of Cambridge, Ph.D. dissertation, p. 61 (2017).
22. Campo, Hannes Marcus Jeff, "Bioengineering strategies of the uterus towards improving current investigative models and female reproductive health". Universitat de València, Ph.D. dissertation, p. 58 (2018).
23. Jarjour, Nicholas Najib, "The Transcription Factor Bhlhe40 Regulates Tissue-Resident Macrophages and Type 2 Immunity". Washington University in St. Louis, Ph.D. dissertation, p. 38 (2019).
24. Legendre, Nicholas, "A Single Allele of MyoD or Myf5 is Necessary for Adult Myogenesis". University of Connecticut Storrs, Ph.D. dissertation, p. 137 (2019).
25. Hodby, Katharine Ailsa, "Investigating the Mechanism of Bone Marrow Failure Observed in Patients with Acute Myeloid Leukaemia". Queen Mary University of London, Ph.D. dissertation, p. 358 (2017).
26. Nguyen, Khiet Tam Christoph, "The role of stromal Hyaluronan in Malignant Melanoma Progression: Investigation in a Has2-knockdown mouse model". Universität Leipzig, Dr. rer. nat. dissertation, p. 44 (2021).
27. Weir, A.F., *Methods Mol. Biol.*, **16**, 1-6 (1993).
28. Weir, A.F., *Methods Mol. Biol.*, **16**, 7-16 (1993).
29. Price, P.A. et al., *J. Biol. Chem.*, **244(4)**, 929-932 (1969).
30. Price, P.A. et al., *J. Biol. Chem.*, **244(4)**, 917-923 (1969).
31. Liao, T.-H., *J. Biol. Chem.*, **250(10)**, 3831-3836 (1975).
32. Lazarides, E., and Lindberg, U., *Proc. Natl. Acad. Sci. USA*, **71(12)**, 4742-4746 (1974).
33. Kunitz, M., *J. Gen. Physiol.*, **33(4)**, 349-362 (1950).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

D4513dat Rev 06/22 VLR,JWM,RBG,GCY,MAM

