

Product Information

# Deoxyribonuclease I from bovine pancreas

Standardized vial containing 2000 Kunitz units of DNase I (D4527), Vial of  $\geq$ 0.25 mg total protein

#### D4263

# **Product Description**

CAS Registry Number: 9003-98-9

Enzyme Commission (EC) Number: 3.1.21.1

Synonyms: DNase I, Deoxyribonuclease A, Deoxyribonucleate 5'-oligonucleotidohydrolase

Deoxyribonuclease I (DNase I) is an endonuclease that cleaves DNA by preferentially acting on phosphodiester bonds adjacent to pyrimidines, to produce polynucleotides with terminal 5'-phosphates. A tetranucleotide is the smallest average digestion product. In the presence of Mg²+ ions, DNase I attacks each strand of DNA independently and the cleavage sites are random. If Mn²+ ions are present, both DNA strands are cleaved at approximately the same site.¹ DNase I hydrolyzes single-stranded DNA, double-stranded DNA, and chromatin (the reaction rate is restricted by DNA association with histones).

DNase I is found in most cells and tissues. In mammals, the pancreas is one of the best sources for the enzyme. Pancreatic DNase I was the first DNase to be isolated. The calculated molecular mass is 30,072 Da. DNase I exists as a mixture of glycoproteins with two disulfide bridges.<sup>2</sup>

Bovine pancreatic DNase I contains four chromatographically distinguishable components, labeled A, B, C, and D.<sup>3</sup> The molar ratios of A:B:C in a pancreatic extract are 4:1:1. Only minor amounts of D are found.

## Carbohydrate Content<sup>4</sup>

Carbohydrate / Form	Α	В	С	
N-Acetylglucosamine	2	3	2	
Mannose	6	5	5	
Sialic Acid	-	1	-	
Galactose	-	1	-	

Forms A and B differ in carbohydrate content (see preceding Table).<sup>4</sup> Form C differs from Forms A and B by having one less His and one more Pro, and in the carbohydrate chain.<sup>4</sup>

DNase I is used to remove DNA from protein and nucleic acid samples, and to nick DNA as a first step to incorporate labeled bases into DNA. Several theses<sup>5-9</sup> and dissertations<sup>10-45</sup> have cited use of product D4263 in their protocols.

## Isoelectric points:2

• A: 5.22

• B: 4.96

• C: 5.06

• D: 4.78

Optimal pH: 7-8

Extinction Coefficient:  $E_{28}\% = 11.1$ 

#### **Activators**

1

- DNase I has an absolute requirement for divalent metal cations.
- The most commonly used divalent metal cation is Mg<sup>2+</sup>.<sup>46,47</sup>
- However, Mn<sup>2+</sup>, Ca<sup>2+</sup>, Co<sup>2+</sup>, and Zn<sup>2+</sup> will activate DNase I.<sup>47-48</sup>
- 5 mM Ca<sup>+2</sup> will stabilize DNase I against proteolytic digestion.<sup>49</sup>
- 0.1 mM Ca<sup>+2</sup> is needed to reduce the rate of inactivation by one-half.<sup>49</sup>



#### **Inhibitors**

There is no general inhibitor specific for DNase I. $^{46,47}$  Citrate inhibits Mg $^{2+}$ -activated DNase I, but not Mn $^{2+}$ -activated DNase I.

- 2-Mercaptoethanol (the reduced enzyme is inactive, but can be reactivated in the presence of Ca<sup>2+</sup> or Mg<sup>2+</sup> ions)<sup>47</sup>
- Chelators (such as EDTA, EGTA)
- Sodium dodecyl sulfate (SDS)<sup>50</sup>
- Actin<sup>51</sup>

## Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

## Product

This product is chromatographically purified from bovine pancreas. The purification procedure is not selective for any form (A, B, C, or D) of DNase I. D4527 is supplied as a lyophilized powder, containing CaCl<sub>2</sub>.

Protein content: ≥ 80% (Biuret)

Specific activity: ≥ 2000 Kunitz units/mg protein

Unit Definition:52

- One Kunitz unit will produce a ΔA<sub>260</sub> of 0.001 per minute per mL at pH 5.0 at 25 °C, using DNA, Type I or III, as substrate, with [Mg<sup>2+</sup>] = 4.2 mM.
- This enzyme assay reaction is performed in 95 mM acetate buffer, pH 5.0, at 25 °C, containing 4.75 mM Mg<sup>2+</sup> and 1.9 mM Ca<sup>2+</sup>, in a 3 mL reaction.

#### Contaminants:

Protease: ≤ 0.005%RNase: ≤ 0.01%

Chymotrypsin: ≤ 0.01%

# Preparation Instructions

This enzyme is soluble in 0.15 M NaCl at 5 mg/mL.

## Storage/Stability

DNase I retains activity for at least three years, when unopened and stored long-term at the recommended temperature, -20 °C.

Solutions of DNase I at 10 mg/mL in 0.15 M NaCl may lose <10% of its activity when stored for a week in aliquots at -20 °C. The same solutions stored in aliquots at 2-8 °C can lose  $\sim20\%$  activity.

DNase I remains active in solution between pH 5-7 up to 60 °C for at least five hours. A 1 mg/mL solution in acetate buffer (pH 5.0) or Tris buffer (pH 7.2) loses activity at the rate of 6% per hour. At 68 °C, DNase I loses activity in <10 minutes.

# References

- Sambrook, J. et al., Molecular Cloning: A Laboratory Manual, 2<sup>nd</sup> edition. Cold Spring Harbor Laboratory Press (Cold Spring Harbor, NY), Vol. 2, p. 5.83 (1989).
- Liao, T.-H. et al., J. Biol. Chem., 248(4), 1489-1495 (1973).
- 3. Kim, H.S., and Liao, T.-H., *Anal. Biochem.*, **119(1)**, 96-101 (1982).
- 4. Salnikow, J. *et al.*, *J. Biol. Chem.*, **245(21)**, 5685-5690 (1970).
- 5. Burton, Jenna Hart, "Regulatory T Cells in Canine Cancer". Colorado State University, M.S. thesis, p. 18 (2011).
- 6. Villanueva, Erika Bianca, "Characterizing the Role of Secretory Phospholipase A<sub>2</sub> Group IIA in Glial Cell-Mediated Neurotoxicity". University of British Columbia, M.Sc. thesis, p. 119 (2011).
- 7. Di Bello, Patrick Louis, "Understanding the Causal Agent of Rose rosette disease". University of Arkansas, M.S. thesis, p. 31 (2015).
- 8. Padmanaban, Veena, "Serum Autoantobodies Block Cancer Invasion in 3D Organotypic Culture". Johns Hopkins University, M.S. thesis, p. 23 (2015).
- 9. Du Plessis, Mar-Dean, "Free-solution Capillary Electrophoresis for the Structural Characterization of Proteins". Western Sydney University, M.Sc. thesis, p. 102 (2018).
- 10. Ellis, Marianne Jane, "Development of a novel hollow fibre membrane for use as a tissue engineered bone graft scaffold". University of Bath, Ph.D. dissertation, p. 52 (2005).
- 11. Heer, Flavio, "CMOS-based Microelectrode Array for Communication with Electrogenic Cells". Swiss Federal Institute of Technology (ETH Zürich), Dr. rer. nat. dissertation, p. 105 (2005).
- 12. Serchov, Tsvetan, "Investigation of photic regulation of the Ras signalling cascade in the visual cortex and circadian systhem of mice". Ruhr University Bochum, Ph.D. dissertation, p. 43 (2007).
- Glick, James J., "DNA Adduction in Context: Native Metabolic Activation of a Foodborne Procarcinogen and Adducted Oligonucleotide Sequencing Monitored by LC-MS/MS". Northeastern University, Ph.D. dissertation, p. 46 (2008).

- 14. Palmer, Lisa Joanne, "Neutrophil extracellular traps in periodontitis". University of Birmingham, Ph.D. dissertation, p. 99 (2010).
- 15. Eisen, Christian, "Development and investigation of a novel model system representing all three subtypes of pancreatic ductal adenocarcinoma reveals novel biomarkers and distinct drug sensitivities". Ruperto-Carola University Heidelberg, Dr. rer. nat. dissertation, p. 37 (2012).
- Gupta, Jalaj, "p38a MAPK in inflammationassociated colorectal cancer". Universidad Autónoma de Madrid, Ph.D. dissertation, p. 67 (2013).
- 17. Illes-Toth, Timea Eva, "Linking the structure of alpha-synuclein oligomers to function in Parkinson's disease". Sheffield Hallam University, Ph.D. dissertation, p. 53 (2013).
- 18. Jorgens, Danielle Mary, "Ultrastructural Characterization of Cellular Interactions Within Resting, Developing and Tumorigenic Breast Epithelium". University of California Berkeley, Ph.D. dissertation, p. 47 (2013).
- 19. Strusi, Valentina, "Toward a 3D *in vitro* model based on decellularized thymus to maintain adult thymic epithelial cells functionality". Università di Bologna, Ph.D. dissertation, p. 91 (2013).
- 20. Kanstinger, Philipp, "Key environmental factors influencing the early life history of clupeoid fishes, adding nutrients and heat". University of Hamburg, Dr. rer. nat. dissertation, p. 65 (2014).
- 21. Markides, Hareklea, "The use of MRI and MNP to image and track cells *in vivo* for arthritic cell-based therapies". Keele University, Ph.D. dissertation, p. 66 (2014).
- Partington, Leanne, "Development of a pressurised transmural decellularisation method for application in tissue engineering trachea". University College London, D.Eng. dissertation, p. 145 (2014).
- 23. Kıray, Hülya, "Identification of Astrocytic Factors That Could Play a Role in Myelination". University of Glasgow, Ph.D. dissertation, p. 47 (2015).
- 24. Lo, Alvin Tu, "Characterization of the Stromal and Epithelial Pathways in the Mammary Gland causing Susceptibility to Cancer". University of California Berkeley, Ph.D. dissertation, p. 23 (2015).
- Nguyen-Ngoc, Kim-Vy, "Novel roles of extracellular matrix and non-muscle myosin ii in regulating epithelial migration and proliferation". Johns Hopkins University, Ph.D. dissertation, p. 56 (2015).

- 26. Scott, Nicole, "Targeting Lysosphingolipid-Induced Lysosomal Impairment as a Critical Point for Therapeutic Intervention in Lysosomal Storage Disorders". University of Rochester, Ph.D. dissertation, pp. 130, 131 (2015).
- 27. Santeramo, Ilaria, "Characterization of Renal CD133+ Cells and their Therapeutic Efficacy in a Model of Acute Kidney Injury". University of Liverpool, Ph.D. dissertation, p. 46 (2016).
- 28. König, Juliana Leonie Berit, "Analyse der Funktionalität von Hormonrezeptoren und der HOXA10-Expression endometrialer Stromazellen zur Ursachenforschung bei Patientinnen mit MRKH-Syndrom" ("Analysis of the functionality of hormone receptors and the HOXA10 expression of endometrial stromal cells for causal research in patients with MRKH syndrome"). Eberhard Karls Universität zu Tübingen, Dr. med. dissertation, p. 104 (2017).
- 29. Zhao, Bing, "The Role of Centrosomal Microtubules in F-actin Dynamics during Neuronal Polarization". University of Hamburg, Ph.D. dissertation, p. 76 (2017).
- Beex, Lana Martina, "Reactive gliosis at the tripartite synapse: Functional characterization of astrocytes in a model of Alzheimer's disease". Universiteit van Amsterdam, Ph.D. dissertation, p. 59 (2018).
- 31. Kong, Joseph Cherng Huei, "Immune Markers to Predict Response to Neoadjuvant Chemoradiotherapy and Identifying Methods to Incorporate Immunotherapy in Locally Advanced Rectal Cancer". University of Melbourne, Ph.D. dissertation, p. 149 (2018).
- 32. Samocha, Alexandr Joseph, "Mammary Epithelial Cell Fate & EGFR-Rasgrp1-Ras Effector Signals". University of California San Francisco, Ph.D. dissertation, pp. 124, 163 (2018).
- 33. Thatenhorst, Denis, "Local Water Fluxes Across the Cell Membrane of Migrating Oligodendrocyte Progenitor Cells". Ruhr University Bochum, Ph.D. dissertation, p. 27 (2018).
- 34. Überbacher, Christa, "Generation of fluorescent reporter lines for hiPSCs derived midbrain dopaminergic neurons: using a CRISPR/Cas9 based knock-in approach to mark TH and GIRK2 expression *in vitro*". Università Degli Studi di Parma, Ph.D. dissertation, p. 161 (2018).
- 35. Bandla, Sravanthi, "Effects of exogenous B-cell immunoglobulin binding protein (BIP) and BIP protein inducer X (BIX) on myelination". National University of Ireland, Galway, Ph.D. dissertation, p. 48 (2019).

- 36. Donella, Hayley Jeton, "Designing Rational Combination Strategies for Overcoming Drug Resistance in Breast Cancer". University of California San Francisco, Ph.D. dissertation, p. 101 (2019).
- 37. Fan, Yuchen, "Vaccine Lipid-Biopolymer Cross-Linked Nanoparticles Against Infectious Diseases and Cancer". University of Michigan, Ph.D. dissertation, p. 94 (2019).
- 38. Prest, Travis Austin, "Development of a Peripheral Nerve Specific Extracellular Matrix-based Hydrogel and its Use in Augmenting Peripheral Nerve Injury and Repair". University of Pittsburgh, Ph.D. dissertation, p. 80 (2019).
- 39. Seelbinder, Benjamin, "Role of Mechanosensitive Chromatin Organization in the Nucleus of Eukaryotic Cells". University of Colorado, Ph.D. dissertation, p. 95 (2019).
- 40. Welle, Anna, "Exploring cell type diversity by comprehensive epigenome and transcriptome analysis". Universität des Saarlandes, Dr. rer. nat. dissertation, p. 35 (2019).
- 41. Fais, Rafael Sobrano, "Papel do NLRP3 na disfunção erétil em modelo de hipertensão DOCA/sal e sua possível ativação por ET-1" ("Role of NLRP3 in erectile dysfunction in DOCA/salt model of hypertension and its possible activation by ET-1"). Universidade de São Paulo, Ph.D. dissertation, p. 61 (2020).
- 42. Saunders, Scott Harrison, "Mechanisms of phenazine-mediated extracellular electron transfer by *Pseudomonas aeruginosa*". California Institute of Technology, Ph.D. dissertation, p. 72 (2020).
- 43. Walker, Sophie, "The Role of iRhom2 in Cardiovascular Pathology". Queen Mary University of London, Ph.D. dissertation, p. 49 (2020).
- 44. Pottast, Maria, "Investigation into the role of the transcription factor RelB for the immunological tolerance induction". Technischen Universität München, Dr. rer. nat. dissertation, p. 31 (2021).
- 45. Wright, Adrienne Beth, "The isolation, culture-expansion, cryopreservation, characterization, and properties of umbilical cord-derived mesenchymal stromal cells and their extracellular vesicles". Kansas State University, Ph.D. dissertation, p. 24 (2021).

- 46. Weir, A.F., Methods Mol. Biol., 16, 1-6 (1993).
- 47. Weir, A.F., Methods Mol. Biol., 16, 7-16 (1993).
- 48. Price, P.A. *et al.*, *J. Biol. Chem.*, **244(4)**, 929-932 (1969).
- 49. Price, P.A. *et al.*, *J. Biol. Chem.*, **244(4)**, 917-923 (1969).
- 50. Liao, T.-H., *J. Biol. Chem.*, **250(10)**, 3831-3836 (1975).
- 51. Lazarides, E., and Lindberg, U., *Proc. Natl. Acad. Sci. USA*, **71(12)**, 4742-4746 (1974).
- 52. Kunitz, M., *J. Gen. Physiol.*, **33(4)**, 349-362 (1950).

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Standard Warranty

The applicable warranty for the products listed in this publication may be found at <a href="SigmaAldrich.com/terms">SigmaAldrich.com/terms</a>.

**Contact Information** 

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

