



Product Information

FIXING SOLUTION 5X Concentrate

Product No. **F 7264**
Store at room temperature

Product Description

This product contains 60% (w/v) trichloroacetic acid and 17.5% (w/v) 5-sulfosalicylic acid. After diluting 5-fold with deionized water, the working solution contains 12% trichloroacetic acid and 3.5% 5-sulfosalicylic acid. It is suitable for fixing proteins in polyacrylamide and agarose gels prior to staining.

Procedure

Protocol for Fixing PAGE, SDS-PAGE and IEF gels

1. Rinse the gel in deionized water for 20-30 seconds to remove the electrode buffer.
2. Place the gel in diluted fixing solution for 30 minutes. Periods longer than 30 minutes are not recommended.
3. Rinse the gel in several changes of deionized water to remove fixing solution completely. Some procedures may require removal of fixing solution by a different method. Consult your particular staining protocol for those requirements.
4. Gel is now ready for staining.

3/99

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.