



Enrichment Sample Homogenizer (ESH)

5623360001

Intended Use:

The Enrichment Sample Homogenizer is used for automatic homogenization of food enrichment samples with enrichment media.

Features of the ESH instrument:

- Multi-function digital display/control panel
- Variable speed (4, 6, 8, and 10 strokes)
- Variable time (1 second to 1 hour or open-running/countdown)
- Close-tight window door
- Adjustable blending power
- Side-by-side paddle stop
- Silent, brushless motor
- Security drip tray
- Removable autoclavable paddles

Category	Specification
Useful volume	50—400 mL
Power source	100V~ to 240V~ 50 to 60 Hz
Size (width x depth x height)	26.5 × 42 × 26 cm 10.433 × 16.45 × 10.24 in.
Weight	16 kg
Blending time	1 second—1 hour, and open running / countdown
Blending speed	4, 6, 8, and 10 strokes



Operator and Equipment Safety

- Place the ESH instrument upright on a clean, flat, stable, horizontal surface, away from any source of excessive heat and close to an easily accessible, properly grounded power supply outlet. Do not run the instrument in any other position
- Do not open the housing without first unplugging the instrument from the power source
- Do not expose the ESH instrument to liquid. If this happens, immediately switch off and disconnect the pump from the power outlet, and then decontaminate the ESH instrument
- Before cleaning, shut down the ESH and disconnect it from the power source
- The power supply must be protected by a fuse below the main connection

For further details, see the Enrichment Sample Homogenizer User Guide

Related Products:

Catalog Number	Instrument
5622900001	DiluCult™ dilutor instrument
5623350001	DiluCult™ 2 dilutor instrument

EDMillipore.com/dilutionpreparation

To place an order or receive technical assistance

In the U.S. and Canada, call toll-free
1-800-645-5476

For other countries across Europe
and the world, please visit:
[EMDMillipore.com/offices](https://www.emdmillipore.com/offices)

For Technical Service, please visit:
[EMDMillipore.com/techservice](https://www.emdmillipore.com/techservice)

[EMDMillipore.com](https://www.emdmillipore.com)

MilliporeSigma
400 Summit Drive
Burlington, MA 01803

