

Quick Reference Guide


PURedit™ Cas9 Protein

PECAS9


Protein Reconstitution

	Recommendations for	
	50 µg vials	250 µg vials
Reconstitute in (µL)	30	50
Concentration (µg/µL)	1.7	5
Concentration (pmol/µL)	10.5	30.8

Upon first use, please read the Product Information Sheet. Actual vial content may vary due to overfill, for exact calculations see the Certificate of Analysis on the product page at SigmaAldrich.com.

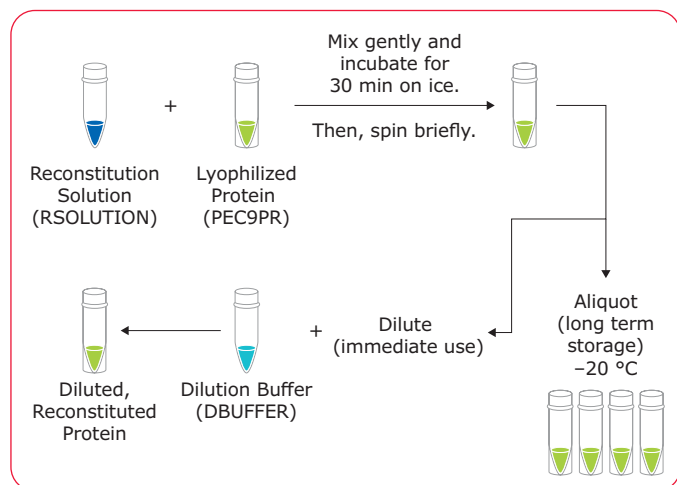

Prepare RNP Complex

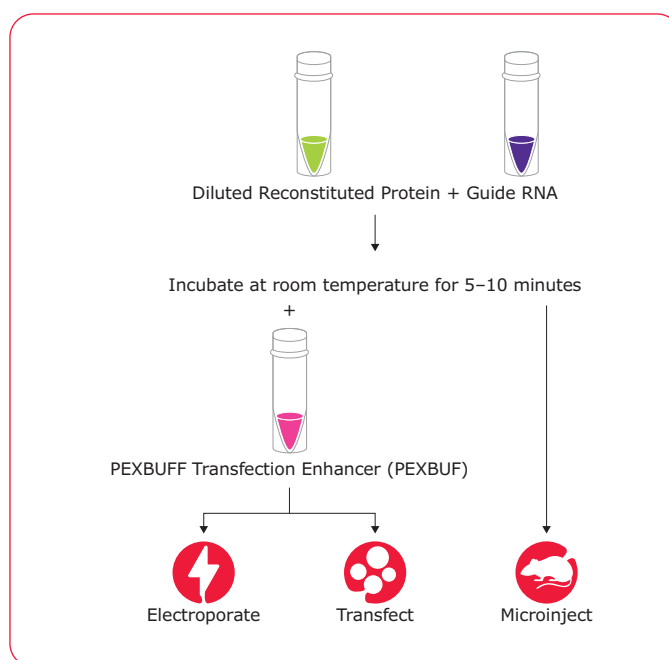

To order guide RNAs, visit
SigmaAldrich.com/CRISPR

Merck, PURedit and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.
 © 2021-2024 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck operates
 as MilliporeSigma in the U.S. and Canada.

MERCK