



Supelco SLB®-IL60 Ionic Liquid GC Columns

Improved Resolution

The variety of analyte-phase interactions possible with the SLB-IL60 column can be leveraged to provide improved separations. This is especially true for complex mixtures comprised of compounds with varying functionality. Polyethylene glycol (PEG) columns rate at 50-52 on our GC Column Polarity Scale; whereas, the SLB-IL60 column has a 60 rating. Therefore, the SLB-IL60 column will exhibit 'wax-like' selectivity for many compound classes. However, the higher rating (60 compared to 50-52) does create subtle selectivity differences which are advantageous for many applications, such as multi-component industrial solvent mixtures.

The SLB-IL60 column was compared directly to five popular commercially available PEG columns, each from a different manufacturer. All columns were 30 m x 0.25 mm I.D., 0.25 μm dimensions, except the SLB-IL60 column, which has a 0.20 μm film thickness. Table 1 shows the maximum temperature limits for all columns tested. Complete specifications of SLB-IL60 columns are shown in Table 2.

Table 1. Maximum Temperature Limits *

Column	Isothermal	Programmed
PEG 1	280 °C	280 °C
PEG 2	260 °C	270 °C
PEG 3	250 °C	260 °C
PEG 4	250 °C	260 °C
PEG 5	280 °C	300 °C
SLB-IL60	300 °C	300 °C

^{*} Obtained from paperwork included with commercial columns.

Table 2. SLB-IL60 Column Specifications

- Application: The SLB-IL60 polar ionic liquid column has a polarity/selectivity similar to that of polyethylene glycol (PEG) columns (usually have 'wax' in the product name), but different enough to provide a unique elution pattern. It also has a higher maximum temperature of 300 °C, compared to 250-280 °C for most PEG columns. These features make it an excellent alternative to existing 'wax' columns. The combination of a high thermal limit and an orthogonal selectivity to non-polar columns also makes it a good GCxGC column choice. Launched in 2012.
- USP Code: None
- Phase: Non-bonded; proprietary
- Temp. Limits: 35 °C to 300 °C (isothermal or programmed)

Industrial Solvents

A 56-component industrial solvent mix that contained alcohols, aldehydes, aromatics, chlorinated hydrocarbons, esters, ethers, ketones, and nitrogen-containing compounds was analyzed on each column under identical conditions. All five PEG columns produced almost identical chromatography. **Figure 1** and **Figure 2** (see back page) show the chromatograms obtained from the PEG 1 and SLB-IL60 columns, respectively. Observations are that the SLB-IL60 column provides:

- An overall quicker analysis (22 minutes compared to 26 minutes)
- A vastly different elution pattern
- Resolution of more analytes; three co-elutions (6 compounds total) compared to six co-elutions (12 compounds total)
- Resolution of a contaminant peak (peak c), suspected to be 2-methylbutyl acetate, from isoamyl acetate (peak 35)

Conclusion

Columns based on polyethylene glycol phase chemistry are widely used for a variety of applications (such as solvents and FAMEs). However, modification of PEG phase chemistry to affect selectivity is very limited. It is advantageous to possess columns with alternative selectivity, because resolution can be greatly affected by selectivity. The SLB-IL60 column is able to undergo many of the same analytephase interactions as PEG columns, plus some additional interactions. This results in the SLB-IL60 column being similar enough to PEG columns to make it useful for many of the same applications, but different enough to impart unique selectivity which can be leveraged to change elution patterns and/or improve resolution.

Featured Products


Description	Cat. No.
SLB-IL60, 15 m x 0.10 mm I.D., 0.08 μm	29503-U**
SLB-IL60, 30 m x 0.25 mm I.D., 0.20 μm	29505-U
SLB-IL60, 60 m x 0.25 mm I.D., 0.20 μm	29506-U**
SLB-IL60, 30 m x 0.32 mm I.D., 0.26 μm	29508-U**
SLB-IL60, 60 m x 0.32 mm I.D., 0.26 μm	29509-U**


^{**}Products will be available soon.

Visit sigma-aldrich.com/il-gc for updates.

For updates, visit sigma-aldrich.com/il-gc



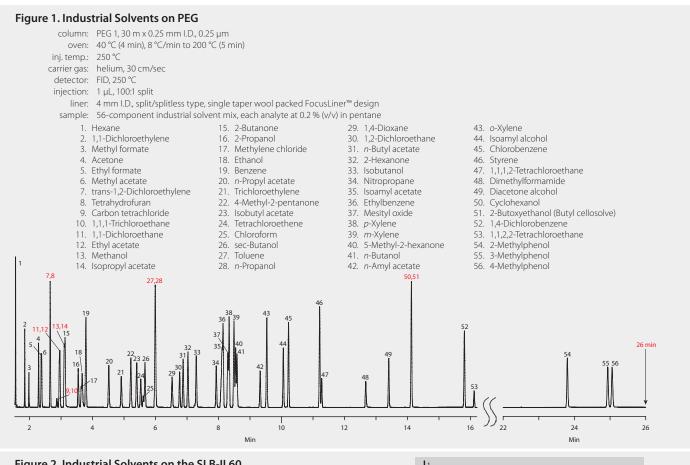

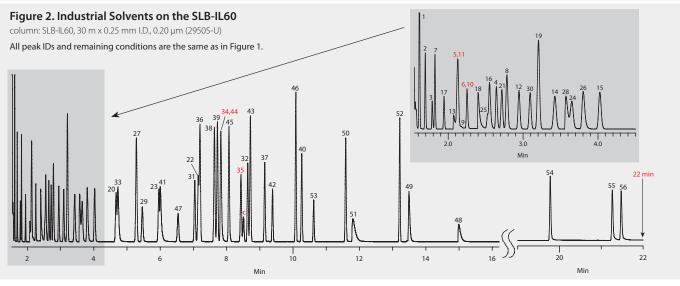

Enabling Science to
Improve the Quality of Life

Order/Customer Service (800) 325-3010 • Fax (800) 325-5052
Technical Service (800) 325-5832 • sigma-aldrich.com/techservice
Development/Custom Manufacturing Inquiries SAFC (800) 244-1173
Safety-related Information sigma-aldrich.com/safetycenter

World Headquarters 3050 Spruce St. St. Louis, MO 63103 (314) 771-5765 sigma-aldrich.com

©2012 Sigma-Aldrich Co. LLC. All rights reserved. SAFC, SLB, SiGMA-ALDRICH and SUPELCO are trademarks of Sigma-Aldrich Co. LLC, registered in the US and other countries. Solutions within is a trademark of Sigma-Aldrich Co. LLC. FocusLiner is a registered trademark of SGE Analytical Science Pty Ltd. Supelco brand products are sold through Sigma-Aldrich, Inc. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see product information on the Sigma-Aldrich website at www.sigmaaldrich.com and/or on the reverse side of the invoice or packing slip.

OSN 11569-509381 / T412091 1092