

User Protocol for 3Y-ZrO₂ Ink for Direct-Ink Writing

Catalog No. [918571](#)

Chemicals Needed:

- 3D Printable Yttria-stabilized Zirconium(IV) Oxide Ink ([918571](#))
- Luperox®, 231, 1,1-Bis(tert-butylperoxy)-3,3,5-trimethylcyclohexane ([479896](#))

Equipment/Consumables Needed:

- Mixer container
- Plastic Spatula
- Balance
- Planetary Mixer
- Syringe (with caps)
- Syringe Nozzle
- Syringe plunger
- Direct-Ink Write (DIW) Printer
- Furnace

Step by Step Procedure for Ink Mixing and Loading into Syringe:

- (1) Place mixer container on balance and tare weight.
- (2) Using the metal spatula, scoop out desired amount of the ink ([918571](#)) from container into the mixer container on the balance. (*For reference, for filling a 10 mL syringe, 30 grams of [918571](#) is sufficient.*)
- (3) Once the desired amount of ink ([918571](#)) is in the mixer container, zero the balance again. Add the Luperox® 231, which will be 0.3 wt% of the total mass of ink ([918571](#)) used. (*For example, if 30 grams of ink ([918571](#)) is used, 0.09 grams of Luperox® 231 will be added.*) A plastic pipette works well to add the Luperox® 231, dropwise into the ink ([918571](#)).
- (4) After the addition of Luperox® 231, use the metal spatula to mix the ink by hand until the paste is visibly uniform, with the viscosity of ink mixture slightly decreasing after the addition.
- (5) Using a planetary mixer, mix the ink using the following mixing sequence:
 - a. Mix setting @ 2000 rpm for 2 min.
 - b. Defoam setting @ 2200 rpm for 3 min.Repeat this mixing sequence twice. With hand mixing in between.
- (6) Once thoroughly mixed, the ink can be loaded into a capped syringe for DIW Printing. Using the metal spatula, scoop the ink from the container and deposit on the side of the syringe. Carefully tap the ink from the syringe wall to fill the syringe fully.

- (7) After all ink is loaded into the syringe. Place the syringe into the mixer and use the Defoam setting @ 2200 rpm for 10 seconds to eliminate trapped air bubbles that will cause inconsistencies in printing.
- (8) After plunger is placed into the syringe, ink is ready for printing.

Other Notes for DIW:

- Syringe Nozzle sizes ranging from 250 μm – 800 μm can be used
- Initial pressure for DIW printing ranges from 12-18 psi, depending on nozzle diameter used

Post Processing of DIW printed parts:

- (1) Curing @ 110°C for 12-16 hours.
- (2) Sintering using the following sequence:
 - Starting at RT, Ramp to 200 °C for 6 hours.
 - Ramp to 300 °C @ 1 °C/min. Hold for 4 hours.
 - Ramp to 400 °C @ 1 °C/min. Hold for 2 hours.
 - Ramp to 800 °C @ 2 °C/min. Hold for 4 hours.
 - Ramp to 1090 °C @ 2 °C/min. Hold for 15 hours.
 - Ramp down to 50 °C @ 2 °C/min.

Copyright © 2021 Merck KGaA, Darmstadt, Germany and/or its affiliates. All rights reserved. Merck, the vibrant M, SigmaAldrich, and TissueFab are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources. More information on our branded products and services on [MerckMillipore.com](https://www.MerckMillipore.com)

The life science business of Merck
operates as MilliporeSigma in the
U.S. and Canada.

MERCK