

**Product Information** 

# SIGMAFAST™ 3,3′-Diaminobenzidine tablets

Tablet, to prepare 5 mL

#### D4293

# **Product Description**

Diaminobenzidine (DAB) is used in many applications for visualization of peroxidase activity. <sup>1-6</sup> DAB is the immunohistology substrate of choice, as it produces an intense brown-black stain, which is resistant to alcohol. Slides stained with DAB may be cover-slipped in the traditional manner and stored for future reference.

SIGMAFAST™ 3,3'-Diaminobenzidine tablets have been developed for use in immunohistology as a precipitating substrate for the detection of peroxidase activity:

Peroxidase + 2  $H_2O_2 \rightarrow O_2 + 2 H_2O$  (pH 7.6)

 $O_2 + DAB \rightarrow insoluble$ , brown-black precipitate

SIGMAFAST™ DAB tablets require no additional ingredients or procedures to prepare an active substrate solution. One SIGMAFAST™ DAB tablet set (one DAB tablet and one Urea Hydrogen Peroxide tablet) dissolved in 5 mL of ultrapure water provides 5 mL of a ready-to-use substrate solution that contains:

- 3,3'-Diaminobenzidine (DAB): 0.7 mg/mL
- Urea Hydrogen Peroxide (H<sub>2</sub>O<sub>2</sub> equivalence, 0.7 mg/mL): 2.0 mg/mL
- Trizma® buffer: 60 mM

This product has been used in studies on such systems as hepatocellular carcinoma, gastric carcinoma, *Xenopus laevis* muscle and embryos, mouse models, 10-14 *Zootoca vivipara* embryos, 5 dogfish testes, 6 rainbow trout, 7 Sprague-Dawley rats, 13 and food samples. 8 Several theses 19-22 and dissertations have cited use of product D4293 in their research protocols.

## Storage/Stability

Store the tablets at -20 °C.

### Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

# Components

- 3,3'-Diaminobenzidine (DAB) tablets (Component Number D9292): 5SET (5 tablets) or 50SET (50 tablets)
- Urea Hydrogen Peroxide tablets (Component Number U1380): 5SET (5 tablets) or 50SET (50 tablets)

# Reagents and Equipment Required but Not Provided

- Ultrapure water (17 MΩ•cm or equivalent)
- Pipette capable of delivering 5 mL
- Test tubes

# Additional Optional Materials

- 0.2 μm filter (such as Cat. No. WHA10462701)
- Nickel(II) chloride (NiCl<sub>2</sub>) or Cobalt(II) chloride (CoCl<sub>2</sub>), to prepare 0.3% (w/v) stock solution to enhance tissue stains
- PBS for washing

1

# **Preparation Instructions**

- 1. Remove the required number of DAB and Urea Hydrogen Peroxide tablets from the freezer.
- 2. Allow the tablets to reach room temperature.
- Open DAB tablet package (silver foil) and Urea Hydrogen Peroxide tablet package (gold foil). Drop the tablets into an appropriate container.
  Do not touch the tablets with your fingers.



- 4. Add 5 mL of ultrapure water.
- 5. Vortex until dissolved.

The SIGMA $FAST^{\text{TM}}$  DAB Substrate Solution is now ready for use. For best results, the solution should be used within one hour.

#### Procedure

- Cover the treated tissue section with 0.2-0.5 mL of DAB Substrate Solution.
- The DAB reaction may occur rapidly. Color development should be carefully monitored during the reaction to prevent overdevelopment and high backgrounds. Reactions may be stopped by gently washing the slide in water or PBS.
- DAB reactions may be enhanced by the addition of a NiCl<sub>2</sub> or CoCl<sub>2</sub> solution. Add 0.5 mL of 0.3% (w/v) stock solution to 4.5 mL DAB Substrate Solution. The addition of metal salts to DAB changes the color of the precipitate product to black or blue-black.
- 4. Occasionally the DAB Substrate Solution may be hazy. The haziness may be removed by filtering the solution through a 0.2 μm filter.
- Tissues stained with SIGMAFAST™ DAB Substrate Solution may be dehydrated with alcohol and mounted with traditional resinous mounting media.

# **Troubleshooting**

### Background is too high

- 1. Use a blocking step prior to the application of the primary antibody. Diluted normal serum (10% v/v) from the same species as the secondary antibody generally produces the best results.
- 2. Block endogenous peroxidase by flooding the slide with a solution of 4 parts methanol and 1 part 3%  $H_2O_2$  solution.
- Decrease the staining time.
- 4. Titer the conjugate to optimize working dilution.

#### No color develops or color is too faint

- 1. Adjust the concentration of the primary antibody.
- Adjust the concentration of the secondary antibody.
- 3. Determine if the enzyme conjugate is active.
- 4. Consider using an amplifying system such as avidin-biotin or peroxidase anti-peroxidase.
- 5. Increase the staining time.

6. Determine if enzymatic treatment (unmasking) of the antigen is required prior to application of the primary antibody.

# References

- 1. Nakane, P.K., and Pierce, G.B., Jr., *J. Histochem. Cytochem.*, **14(12)**, 929-931 (1966).
- Trojanowski, J.Q. et al., J. Histochem. Cytochem., 31, 1217-1223 (1983).
- DeJong, A.S.H. et al., Histochem. J., 17(10), 1119-1130 (1985).
- Chu, N.M. et al., J. Histochem. Cytochem., 37(2), 257-263 (1989).
- Merchenthaler, I. et al., "Silver Intensification in Immunocytochemistry", in Techniques in Immunocytochemistry (Bullock, G., and Petrusz, P., eds.). Academic Press Ltd. (San Diego, CA), pp. 217-252 (1989).
- 6. Hsu, S., and Soban, E., *J. Histochem. Cytochem.*, **30(10)**, 1079-1082 (1982).
- Yarchoan, M. et al., Clin. Cancer Res., 23(23), 7333-7339 (2017).
- 8. Tsamandas, A.C. et al., Anticancer Res., **29(2)**, 703-709 (2009).
- Lunardi, A., and Dente, L., *Mech. Dev.*, **119** Suppl 1, S49-S54 (2002).
- 10. Nolan-Stevaux, O. et al., Genes Dev., **23(1)**, 24-36 (2009).
- 11. Brotherton, T.E. *et al.*, *Proc. Nat. Acad. Sci. USA*, **109(14)**, 5505-5510 (2012).
- van Dop, W.A. et al., Gastroenterology, 136(7), 2195-2203 (2009).
- 13. Nandi, S.S. *et al.*, *Diabetes*, **65(10)**, 3075-3090 (2016).
- 14. Wakabayashi, T. *et al.*, *Arterioscler. Thromb. Vasc. Biol.*, **38(11)**, 2576-2589 (2018).
- 15. Stewart, J.R. *et al.*, *J. Exp. Biol.*, **214(18)**, 2999-3004 (2011).
- 16. Gautier, A. et al., *Biol. Reprod.*, **91(4)**, 1-15 (2014).
- 17. Jørgensen, T.R. *et al.*, *Folia Parasitol.* (*Praha*), **56(4)**, 251-258 (2009).
- 18. Boutten, B. *et al.*, *Food Agr. Immunol.*, **11(1)**, 51-59 (1999).
- 19. Bhatla, Sunetra, "Quantification and RNAi knockdown phenotype of *RNP-4F* isoforms during CNS development in *Drosophila*". Miami University, M.S. thesis, p. 72 (2006).
- Goolia, Melissa Nalini, "The Expression and Localization of the V and W Accessory Proteins in Porcine and Human Cells Infected with Nipah Virus". University of Manitoba, M.Sc. thesis, pp. 39, 40, 42, 43 (2009).



- 21. Coutinho, Frazer Paul, "An investigation of CD147 expression in apoptotic vesicles". University of Otago, M.Sc. thesis, p. 20 (2013).
- 22. Springer, Isaac C., "Investigation into the Role of Platelet Derived Growth Factor (PDGF) in Type I Collagen Glomerulopathy". Missouri State University, M.S. thesis, p. 38 (2018).
- 23. Balaskas, Christos, "Cellular development of the enteric nervous system in the chick embryo". University College London, Ph.D. dissertation, pp. 50, 73, 222 (1997).
- 24. McFetridge, Peter Stuart, "The use of porcine carotid arteries as a matrix material for tissue engineered small diameter vascular grafts". University of Bath, Ph.D. dissertation, p. 53 (2002).
- 25. Bungay, Anton, "T<sub>3</sub> as a primary hepatic mitogen". University College London, Ph.D. dissertation, p. 52 (2008).
- 26. Eozenou, Caroline, "FOXL2: A regulator of endometrial physiology? First insights from ruminants". Université Paris-Sud, Ph.D. dissertation, pp. 192, 220 (2013).
- 27. Siew, Leonard Quok Chean, "The 'smoking endotype' of asthma: studies on cigarette smoke-induced, IL-17A dependent neutrophilic inflammation in the bronchial mucosa of asthmatics who smoke". King's College London, Ph.D. dissertation, p. 69 (2014).
- 28. Filatova, Anna, "The role of Fam20a in the generation of Amelogenesis Imperfecta". Universität Zürich, Dr. sc. nat. dissertation, p. 33 (2015).
- 29. Du, Shengnan, "Novel Microtubule-Disrupting Indole-Based Chalcones That Induce Cell Death in Glioblastoma". University of Toledo, Ph.D. dissertation, p. 110 (2017).
- 30. Erbeldinger, Nadine, "Microvascular Damage as Initial Event of Scar Formation after Carbon Ion Irradiation of Cardiac Substructures". Technischen Universität Darmstadt, Dr. rer. nat. dissertation, p. 26 (2017).
- 31. Urakova, Nadezda, "Identifying molecular virulence factors of rabbit haemorrhagic disease virus". University of Canberra, Ph.D. dissertation, p. 15 (2017).
- 32. Sheikh, Imran S., "The Role of C3-C4 Propriospinal Interneurons on Reaching and Grasping Behaviors Pre- and Post-Cervical Spinal Cord Injury". Temple University, Ph.D. dissertation, pp. 66, 108 (2018).

- Hussain, Ainy Khan, "Non-Immune Islet Beta Cell Susceptibility to Failure in NOD<sup>K</sup> Mice". Australian National University, Ph.D. dissertation, p. 67 (2017).
- 34. Kresinsky, Anne, "Receptor protein-tyrosine phosphatases controlling activity of the oncoprotein FLT3 ITD". Friedrich-Schiller-Universität Jena, Dr. rer. nat. dissertation, p. 45 (2019).
- 35. Syafiq, Tengku Muhamad Faris, "Modeling human genetic kidney malformations caused by *HNF1B* mutation using pluripotent stem cells". University of Manchester, Ph.D. dissertation, p. 63 (2019).
- 36. Reolizo, Lien Mari P., "Potential of Proline Rich Homeodomain (PRH) to inhibit intimal thickening". University of Bristol, Ph.D. dissertation, pp. 53-56 (2021).

### **Notice**

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

#### Technical Assistance

Visit the tech service page at <u>SigmaAldrich.com/techservice</u>.

#### Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

#### **Contact Information**

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

