

Protocol

TissueFab® bioink kit, (Gel)ma Laminin – Vis/405 nm, low endotoxin

Protocol for Catalog No. [926000](#)

Introduction

The TissueFab® bioink kit (Gel)ma Laminin Vis/405 nm features a ready-to-use bioink formulated to delivery exceptional cell viability and printability. Specifically designed for extrusion-based 3D bioprinting, this ready-to-use bioink is further enhanced through 405 nm light crosslinking and the incorporation of laminin, optimizing its performance for advanced tissue fabrication. Compatible with most extrusion-based bioprinters, these biodegradable bioinks are well-suited for use with human mesenchymal stem cells (hMSCs) and various other cell types. The TissueFab® bioink kit (Gel)ma Laminin Vis/405 nm enables the precise creation of 3D cell models and tissue constructs, offering unparalleled potential for research in 3D cell biology, tissue engineering, in vitro tissue models, and regenerative medicine.

Disclaimer

TissueFab® bioink kit (Gel)ma Laminin Vis/405 nm is for research use only; not suitable for human, animal, or other use. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Specifications

Storage	Store TissueFab® bioink kit (Gel)ma Laminin Vis/405 nm at -20 °C . Protect from light by storing the bottle in a foil bag or wrapping it in aluminum foil.
Stability	Refer to the expiration date on the batch-specific Certificate of Analysis.

Materials

Materials supplied

TissueFab® - GelMA-Laminin bioink kit Vis/405 is supplied as follows:

Catalog Number	Quantity
918741	1 x 10ml (1 unit)
L2020	1 x 1ml (1 unit)

Materials required, but not supplied

- Cultured cells (visit our website for an up-to-date list of cell types) link: <https://www.sigmaaldrich.com/life-science/cell-culture/mammalian-cell-lines.html>
- Appropriate cell culture medium
- PBS (Cat. No. [D8537](#))
- Sterile pipette tips for transferring bioink
- Sterile printing cartridge, piston, and nozzle/needle for 3D printing
- Extrusion-based 3D bioprinter
- Water bath or incubator
- Micropipettes
- 405 nm light source



Before you start: Important tips for optimal bioprinting results

Optimize printing conditions. Optimize printing conditions (e.g., nozzle diameter, printing speed, printing pressure, temperature, cell density) for the features of your 3D printer and for your application to ensure successful bioprinting. The suggestions below can guide you.

Reduce bubble formation. If the bioink has air bubbles, the bubbles may hamper bioprinting. Carefully handle the bioink when you mix and transfer it to avoid bubble formation. Do not vortex or shake vigorously.

Aseptic techniques. Follow standard aseptic handling techniques when you prepare and print the bioink, and during cell culture.

Cell density. Resuspend the cell pellet to the appropriate volume for the desired printed structure and cell density. Typical cell density for extrusion-based bioprinting is 1 to 5×10^6 cells/mL. For example, Human bone marrow derived mesenchymal stem cells (hMSCs) have been printed with TissueFab® - GelMA Laminin bioink kit Vis/405 at a concentration of 5×10^6 cells/mL.

Note: The number of prints obtained from each 10-mL bottle of bioink (a unit) will vary depending on the structure that is printed. For example, each 10-mL bottle contains enough material to print a 30- μ L structure in each well of three 96-well plates or a 100- μ L structure in each well of four 24-well plates.

Procedure

A. Prepare bioink kit components

1. Warm the bottle of TissueFab® - (Gel)ma bioink Vis/405 nm in a water bath or incubator set to 37 °C for 30-60 minutes or until the bioink becomes fluid so that it is easy to pipette.
2. Warm the bottle of Laminin in a water bath or incubator set to 37 °C for 5-15 minutes or until the solution has become fluid.

B. Prepare bioink-cell solution

1. centrifuge the cell suspension to obtain a cell pellet. Remove the supernatant carefully so that the cell pellet is not disrupted. Excess supernatant will dilute the bioink and may limit printability.
2. Resuspend the cell pellet at the desired cell density with the bioink kit components according to the table below:

Final Laminin Concentration (μ g/mL)	Bioink Volume (mL)	Laminin Volume (mL)	Media Volume (mL)
0	0.8	0	0.2
50	0.8	0.05	0.15
100	0.8	0.1	0.1
200	0.8	0.2	0.2

3. Gently and slowly pipette the bioink-cell solution up and down several times to ensure the cells are evenly distributed in the bioink solution. Avoid creating air bubbles. DO NOT vortex or shake vigorously.
4. Pipette the bioink-cell solution into the desired printing cartridge. This step creates a filled printing cartridge.
5. Place the remaining bioink in a foil bag or wrap in aluminum foil and store at -20 °C to protect from heat and light.

C. Bioprint

1. Cool the filled printing cartridge to 15-20 °C using a “temperature-controlled printhead”, if available, or place the cartridge in a 4 °C refrigerator for 10–15 minutes to induce gelation.
2. Follow the manufacturer’s 3D printer instructions. Load the print cartridge onto the 3D printer and print directly onto a Petri dish or into multi-well plates. Adjust the flow rate according to the nozzle diameter, printing speed, printing pressure, and temperature.



Example

Printer: Cellink BIO X™ or Cellink INKREDIBLE™ printer

Temperature: 20 °C

Flow rate (speed): 10 mm/s

Nozzle: 22G TT tapered needle

Pressure: 70-90 kPa

D. Crosslink

Place the light source directly above the 3D-bioprinted structure and expose the structure to light (recommended settings: wavelength – 405 nm; irradiance – 10 mW/cm²; exposure – 30-60s). Use the appropriate distance and exposure time based on your light source. For low-intensity light sources, usually available in desktop bioprinters, such as Cellink™ bioprinters (Bio X™ and INKREDIBLE™ printers), distances of 3 cm or less and exposure times of 60s or more are recommended.

The 3D-bioprinted structure is ready for culture or analysis immediately after crosslinking is done.

E. Culture cells.

Culture the bioprinted tissue with the appropriate cell culture medium following standard tissue culture procedures.

Troubleshooting

1. Bioink is incubated at 37°C for 30 minutes, but it is still gel.

Possible reasons – Malfunction of the incubator; bioink is crosslinked due to light exposure.

Solution – Make sure the temperature of the incubator/water bath is correct and make sure the bioink bottle is properly and evenly heated in the incubator/water bath. Do not expose the bioink to light before printing.

2. Air bubble is trapped in the middle of the bioink in the cartridge.

Possible reason – Air bubble was created during transferred or when cells were dispersed in the bioink.

Solution - Warm the cartridge at 37°C for 5–10 minutes or until the bioink becomes fluid. Turn the cartridge so that the tip faces up to allow any air bubbles to exit from the tip of the cartridge. Gently tap the cartridge to help the air bubbles pass through the tip.

3. Printed structure spreads and does not hold its shape.

Possible reasons – Bioink was diluted with cell culture medium that remained in the cell pellet; bioink was not cooled sufficiently before printing, or the printing pressure is too high.

Solution – Do not dilute the bioink. Make sure the bioink has been cooled according to the instructions before printing. Adjust printing pressure to achieve sufficient flow of bioink.

4. Interrupted flow or no flow during printing.

Possible reason – Insufficient printing pressure or nozzle is partially or fully clogged.

Solution – Adjust the printing pressure to achieve sufficient flow of bioink. If the problem persists, change the nozzle.

5. Printed structure dissolves in cell culture medium.

Possible reason – Insufficient crosslinking; exposure to incorrect wavelength; malfunction of the light source.

Solution – Make sure that the light source has sufficient power output and that the printed structure is exposed to the correct wavelength for the appropriate exposure according to the instructions.



Related Products

Name	Cat. No.
TissueFab® - GelAlg-UV bioink	905410
TissueFab® - GelAlg-Vis bioink	906913
TissueFab® - GelMA-UV bioink	905429
TissueFab® - Sacrificial bioink	906905
TissueFab® - Bone Support bioink	915637
TissueFab® - GelMA-Bone-UV bioink	915025
TissueFab® - GelMA-Bone-Vis bioink	915033
TissueFab® - GelMA-Conductive-UV bioink	915726
TissueFab® - GelMA-Conductive-Vis bioink	915963
TissueFab® - Crosslinking solution	919926
TissueFab® - HAMA-UV bioink	919632
TissueFab® - HAMA-Vis bioink	919624
TissueFab® - GelMA-Vis-LAP bioink, low endotoxin	918741
TissueFab® - AlgMA-UV bioink	920983
TissueFab® - AlgMA-Vis bioink	921610
TissueFab® - HAMA-AlgMA-UV bioink	920975
TissueFab® - HAMA-AlgMA-Vis bioink	922862

