

User Manual

ES Cell Characterization Kit

For 100 Tests

SCR001**FOR RESEARCH USE ONLY****Not for use in diagnostic procedures. Not for Human or Animal Consumption.**

Product Overview

Stem cells have become the subject of extensive investigation recently, partly due to their therapeutic potential and because they raise several fundamental issues concerning the regulation of proliferation and differentiation.

Embryonic stem (ES) cells are derived from totipotent cells of the early mammalian embryo and are capable of unlimited, undifferentiated proliferation in vitro.^{1,2,3} Undifferentiated mouse ES cells can be maintained for a long time in media containing the cytokine, leukemia-inhibitory factor (LIF) or our proprietary ES cell culture reagent, ESGRO®.^{4,5} However, upon removal of LIF from the culture medium, in vitro, the mouse ES cells start to differentiate into cells derived from all three germ layers. In contrast, human ES cultures require mouse fibroblasts as feeder cells and cannot be maintained with LIF for self-renewal.^{6,7}

The undifferentiated state of the embryonic stem cell is characterized by high level of expression of alkaline phosphatase (AP)⁸ and the stem cell transcription factor, Oct-4. Nevertheless, these ES cells also exhibit marked differences from their murine counterparts in regard to their expression of stage-specific embryonic antigen (SSEA), that typify undifferentiated human ES and embryonic carcinoma (EC) cells.

SSEA-1, a carbohydrate antigen, is a fucosylated derivative of type 2 polylactosamine and appears during late cleavage stages of mouse embryos. It is strongly expressed by undifferentiated, murine ES cells.^{9,10} Upon differentiation, murine ES cells are characterized by the loss of SSEA-1 expression and may be accompanied, in some instances, by the appearance of SSEA-3 and SSEA-4.¹¹ In contrast, human ES and EC cells typically express SSEA-3 and SSEA-4 but not SSEA-1, while their differentiation is characterized by downregulation of SSEA-3 and SSEA-4 and an upregulation of SSEA-1.^{12,13} Undifferentiated, human ES cells also express the keratin sulphate-associated antigens, TRA-1-60 and TRA-1-81.¹⁴

Our ES Cell Characterization Kit (SCR001) is a specific and sensitive tool for the phenotypic assessment of the differentiation status of ES cells by measuring their AP activity, cell-surface stage-specific antigens (SSEA-1, SSEA-4) as well as expression of TRA-1-60, TRA-1-81 antigens.

Antibody Specificity

	SSEA-1	SSEA-4	TRA-1-60	TRA-1-81
Human ES cell	×	✓	✓	✓
Murine ES cell	✓	×	×	×

Materials Provided

- Fast Red Violet solution (0.8 g/L stock) (90239): Two 15 mL bottles
- Naphthol AS-BI phosphate solution (2 mg/mL) in a buffer, pH 8.5 (CS235583): Two 15 mL bottles
- MS X SSEA-1, IgM, clone MC-480 (90230): One vial containing 100 µL of 1 mg/mL monoclonal antibody
- MS X SSEA-4, IgG, clone MC-813-70 (90231): One vial containing 100 µL of 1 mg/mL monoclonal antibody
- MS X TRA-1-60, IgM, clone TRA-1-60 (90232): One vial containing 100 µL of 1 mg/mL monoclonal antibody
- MS X TRA-1-81, IgM, clone TRA-1-81 (90233). One vial containing 100 µL of 1 mg/mL monoclonal antibody

Materials Required (Not supplied)

- Fixative (e.g., 4% Paraformaldehyde)
- 1X Rinse Buffer (e.g., TBST: 20 mM Tris-HCl, pH 7.4, 0.15 M NaCl, 0.05% Tween®-20)
- Blocking Solution
- Fluorescent-labeled secondary antibodies
- Hematoxylin
- Microscope

Storage and Stability

The ES Cell Characterization Kit consists of two components used for alkaline phosphatase activity determination as well as four ES cell-specific antibodies required to perform 100 tests (including controls).

When stored at 2 to 8 °C, the kit components are stable up to the expiration date. Do not freeze or expose to elevated temperatures. Discard any remaining reagents after the expiration date.

Protocol

Preparation of Reagents

Naphthol/Fast Red Violet Solution

Mix Fast Red Violet (FRV) with Naphthol AS-BI phosphate solution (2 mg/mL) in a buffer, pH 8.5 in a 1:1 ratio (FRV:Naphthol).

Staining Protocol

Alkaline Phosphatase Staining Procedure

1. Culture ES cells for five days prior to analyzing AP activity, at low to medium density.
Note: This time-period is critical if activity levels of AP needs to be observed. According to our findings, five days of culturing are optimal for good AP stain visualization.
2. On day five, aspirate media and fix the ES or EC cells with a fixative (e.g., 4% paraformaldehyde in PBS) for 1-2 minutes.
Note: Do not overfix. Fixing cells longer than 2 minutes will result in the inactivation of alkaline phosphatase.
3. Aspirate fixative and rinse with 1X Rinse Buffer. DO NOT allow wells to dry.
4. Prepare reagents for Alkaline Phosphatase staining as described in [Preparation of Reagents](#) section.
5. Add enough stain solution to cover each well (e.g., 0.5 mL for a well of a 24-well plate). Incubate in dark at room temperature for 15 minutes.
6. Aspirate staining solution and rinse wells with 1X Rinse Buffer. Cover cells with 1X PBS to prevent drying and then count the number of colonies expressing AP (red stem cell colonies), versus the number of differentiated colonies (colorless).
AP staining criteria: Greater than 90% of colonies should remain undifferentiated and express alkaline phosphatase in the well containing 10^3 Units of LIF. P value shall be ≥ 0.05 .

Immunofluorescence Staining Procedure

The cell-surface-antigen expression of cultured cells can be analyzed by using immunofluorescence techniques. The following primary monoclonal antibodies are used to detect surface-antigen expression: anti-SSEA-1, anti-SSEA-4, TRA-1-60 and TRA-1-81 (provided in kit). Fluorescein isothiocyanate (FITC)-labeled goat anti-mouse IgM or anti-IgG can be used as secondary antibody, appropriate to the isotype of the primary antibody.

1. Fix cultured ES cells in 4% paraformaldehyde/PBS for 15-20 minutes at room temperature.
2. Wash twice (5-10 minutes each) with 1X Rinse Buffer.
3. Permeabilize cells with 0.1% Triton® X-100/PBS for 10 minutes at room temperature.
4. Wash twice (5-10 minutes each) with 1X Rinse Buffer.
5. Apply a Blocking solution (e.g., 4% normal goat serum/PBS) for 30 minutes at room temperature.
6. Dilute primary antibodies to working concentrations, in blocking solution (SSEA-1, SSEA-4, TRA-1-60 and TRA-1-81 can be used in the range of 1:10 – 1:50). Incubate primary antibodies for 1 hour at room temperature.
7. Wash three times (5-10 minutes each) with 1X Rinse Buffer.
8. Dilute secondary antibodies in 1X PBS just before use. Incubate secondary antibodies for 30-60 minutes at room temperature.
9. Wash three times (5-10 minutes each) with 1X Rinse Buffer.
10. If stained in plate wells, cells should be covered with 1X PBS for visualization. However, if cells are stained on a coverslip, mount on a slide by using antifade mounting solution.
11. Fluorescence images can be visualized with a fluorescence microscope.
Note: Be sure to use the correct filter when visualizing fluorescently labeled cells.

Product Performance

Staining results obtained with ES Cell Characterization Kit

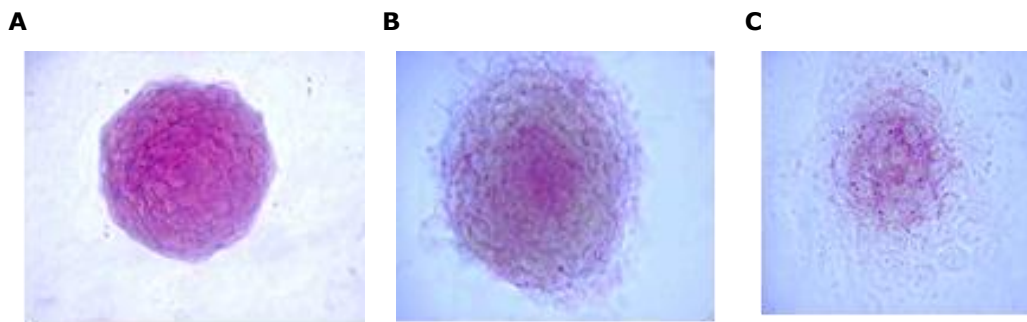


Figure 1: Alkaline Phosphatase staining of ES cells. High magnification revealed (A) Undifferentiated ES cells (mouse MBL.5 cell line) – cultured for five days in media containing LIF/ESGRO®. A concentration of 10^3 units/mL is used for inhibition of differentiation. (B) Differentiated ES cells – cultured at low-medium density for three days in media without any LIF/ESGRO®. (C) Differentiated ES cells – cultured at low-medium density for six days in media without any LIF/ESGRO®.

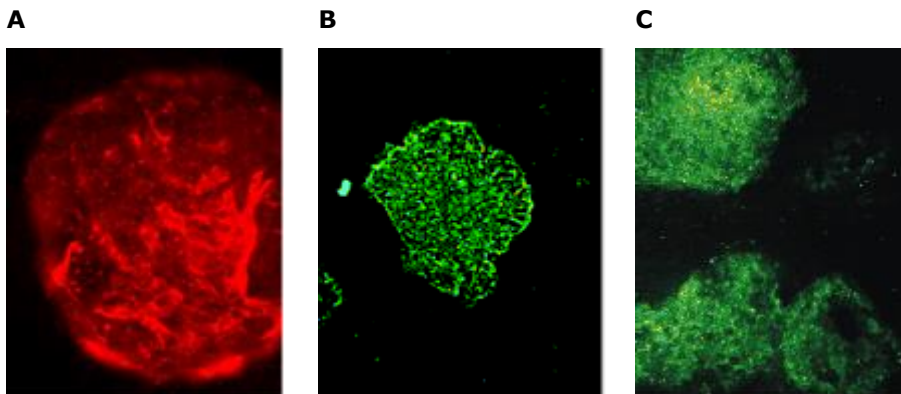


Figure 2: Immunofluorescence images of ES cells, stained with anti-SSEA and anti-TRA antibodies. (A) Undifferentiated, murine ES cells (cultured in presence of LIF) revealed positive immunostaining for SSEA-1. (B) Immunofluorescence staining of human ES cell colony with anti-SSEA-4 monoclonal antibody. As anticipated, ES cells within the colony exhibit strong immunoreactivity to this antibody. (C) Immunofluorescence staining of human ES cells with anti-TRA-1-60 monoclonal antibody. As revealed by the image, the ES cells within the colony exhibit strong immunoreactivity to this antibody.

Photographs B and C courtesy of Dr. Jonathan Draper, University of Sheffield, Department of Biomedical Science.

Related Products

The following related products are available from [SigmaAldrich.com](https://www.sigmaaldrich.com) as separate items.

- SSEA-1 Monoclonal Antibody, purified 100 µg (MAB4301)
- SSEA-3 Monoclonal Antibody, purified 100 µg (MAB4303)
- SSEA-4 Monoclonal Antibody, purified 100 µg (MAB4304)
- TRA-1-60 Monoclonal Antibody, purified 100 µg (MAB4360)
- TRA-1-81 Monoclonal Antibody, purified 100 µg (MAB4381)
- Murine LIF, 5 µg (LIF2005)
- Murine LIF, 10 µg (LIF2010)
- ESGRO®, 1 x 10⁶ units (ESG1106)
- ESGRO®, 1 x 10⁷ units (ESG1107)

References

1. Evans M, Kaufman M. Nature 292, 154 (1981).
2. Martin G. Proc Natl Acad Sci USA 78, 7634 (1981).
3. Morrison SJ, Shah NM, Anderson DJ. Cell 88, 287-298 (1997).
4. Smith AG, Heath JK et al. Nature 336, 688-690 (1988).
5. Williams RL, Hilton DJ et al. Nature 366, 684-687 (1988).
6. Thomson JA, Marshall VS. Curr Top Dev Biol 38, 133 (1998).
7. Shambloott MJ, Axelman J et al. Proc Natl Acad Sci USA 95, 13726-13731 (1998).
8. Pease S, Braghetta P et al. Develop Biol 141, 344 (1990).
9. Solter D, Knowles BB. Proc Natl Acad Sci USA 75, 5565-5569 (1978).
10. Gooi HC, Feizi T, Kapadia A et al. Nature 292, 156-158 (1981).
11. Solter D, Shevinsky LH, Knowles BB. Dev Biol 70, 515-521 (1979).
12. Andrews PW. Dev Biol 103, 285-293 (1984).
13. Fenderson BA, Andrews PW et al. Dev Biol 122, 21-34 (1987).
14. Andrews PW, Banting GS et al. Hybridoma 3, 347-361 (1984).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates
as MilliporeSigma in the U.S. and Canada.

Merck, ESGRO and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2009-2025 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

Document Template 20769660 Ver 4.0

SCR001 Ver 6.0, Rev 20MAY2025, CJ

