Safety Sheet

Heater Shaker Module

Select language:

English	2	Italiano (Italian)	30
العربية (Arabic)	4	日本語 (Japanese)	32
Български (Bulgarian)	6	한국어 (Korean)	34
中文 (Chinese, Simplified)	8	Latviski (Latvian)	36
Hrvatski (Croatian)	10	Lietuvių (Lithuanian)	38
Čeština (Czech)	12	Malti (Maltese)	40
Dansk (Danish)	14	Polski (Polish)	42
Nederlands (Dutch)	16	Português (Portuguese)	44
Eesti keel (Estonian)	18	Română (Romanian)	46
Suomi (Finnish)	20	Srpski (Serbian)	48
Français (French)	22	Slovenčina (Slovak)	50
Deutsch (German)	24	Slovenščina (Slovenian)	52
Ελληνικά (Greek)	26	Español (Spanish)	54
Magyar (Hungarian)	28	Svenska (Swedish)	56

Intended Use

The Opentrons Heater-Shaker Module GEN1 provides on-deck heating and orbital shaking. The Heater-Shaker can heat samples to 95 °C and shake them at speeds ranging from 200 to 3000 rpm. It is compatible with the Opentrons Flex and OT-2 liquid handling robots and selected flat, deep-well, and 96-well plates. The Heater-Shaker can also be used alongside other Opentrons modules and with the Opentrons Flex Gripper.

Specifications

Power Requirements

Power Adapter

Input voltage	100-240 VAC (+/- 10%)
Input frequency	50/60 Hz
Input current	4 A/115 VAC 2 A/230 VAC
Output voltage	36 VDC
Output current	6.1 A
Output wattage	219.6 W maximum
Power source Required	C13

Instrument

Outer dimensions	Length	152 mm (5.98 inches)
	Width	90 mm (3.54 Inches)
	Height	82 mm (3.2 inches)
Weight		1.34 kg

Environmental Conditions

For indoor use only. This module is not rated for wet locations or any other outdoor parameter. See the manual for recommended parameters for system operation.

	Acceptable for Operation	Storage/Transport
Ambient Temperature	20 °C to 24 °C 68 °F to 75 °F	-10 °C to 60 °C 14 °F to 140 °F
Relative Humidity	30-80% non-condensing (below 30 °C)	10-85% non-condensing (below 30 °C)
Altitude	Up to 500 m above sea level	Up to 2000 m above sea level

IEC61010-1 Pollution Degree II

Overvoltage Category II

Ingress Protection Rating IP2X

⚠ Precautions

Please read these safety instructions and manual carefully before using this module. Keep this Safety Sheet near the instrument for reference.

- This module generates enough heat to cause serious burns. Allow maximum clearance to avoid burns. Do not touch if hot.
- If the Status Light is red or amber, avoid touching the module. A white light indicates the module may be cool enough to touch.
- If the power is off, wait at least 10 minutes from last use before touching.
- Safety goggles should be worn at all times during operation.
- Unplug when not in use.

Electrical Warnings

Failure to follow any of these warnings may result in injury, fire, and electric shock.

- Only use the power cord provided.
- Do not use a power cord with inadequate current or voltage ratings. Do not replace the power cord unless at the direction of Technical Service.
- Keep the power cord free of obstructions so you can remove it if necessary.
- Do not connect (plug in), disconnect (unplug), or use AC power cables if:
 - The cable is frayed or damaged.
 - Other attached cables, cords, or receptacles are frayed or damaged.
- Using damaged power cords can cause an electric shock hazard resulting in serious injury or damage to the robot.
- Do not replace the power cable unless at the direction of Opentrons Support.
- Do not attempt to repair or remove the outer case of the module or its power supply unless directed by Opentrons Support.

Maintenance

- Power OFF and disconnect any USB or power cables before beginning cleaning.
- No internal parts require cleaning. The instrument should never be disassembled for cleaning.
- Gently wipe the module surfaces with a solution of 70% diluted alcohol (ethyl/ethanol, isopropyl or methanol) and distilled water. Do not use 100% alcohol. Air dry.
- Do not use acetone. This may cause damage or dissolve parts of the instrument.
- Do not autoclave.
- See the manual Cleaning section for detailed cleaning instruction.

Technical Service

Do not attempt to modify, revise, or repair the instrument. Only authorized Opentrons Support may service the instrument. Unauthorized repair or service may void the warranty and/or cause hazardous conditions.

Installation

- Read the manual for detailed installation directions before installing any module.
- Ensure the module is properly oriented in the caddy sized for this module.
- Tighten the caddy anchors. The shaking motion requires anchoring to prevent the module from dislodging itself while in operation.
- Firmly connect all cords to the module, as directed in the manual.
- The mains power cord and plug must be easily accessible in the case of a hazard in which the cord and plug need to be pulled to remove power from the product.

Equipment Operation

Read the manual before operating.

Power ON/OFF

Plug module into the wall outlet. From the front of the caddy, press (-) on the left side of the power switch for ON. A light (left of the switch) will turn on.

To turn off, press (o) on the right side of the power switch. The light will turn off. Unplug power cord from outlet when not in use.

Emergency Disconnect

Push down the Emergency Stop Pendant (E-stop) to quickly stop robot motion. The robot requires a functional, disengaged E-stop to be attached at all times. The E-stop does not cut power.

In case of an emergency, unplug the power cord from the power source. Contact Opentrons Support before restarting the robot.

Symbol Definitions

	CAUTION, documentation must be consulted
	CAUTION, possibility of electrical shock
	CAUTION, hot surface
	CAUTION, risk of pinching
	CAUTION, risk of explosion
	Manufacturer/ Legal entity
	This product has been evaluated by an OSHA certified Nationally Recognized Testing laboratory (NRTL) in the United States
	Conforms to UL Std 61010-1
	Certified to CSA Std 61010-1
	European Council Directives conformity marking
	United Kingdom conformity marking

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© OPENTRONS 2025
Opentrons Labworks, Inc.
Opentrons Flex™ (Opentrons Labworks, Inc.) Registered names, trademarks,